

Instructions for Use

APP betaCTF ELISA

Enzyme immunoassay for the quantitative determination of human APP β CTF in cell lysate.

REF **JP27776**

 **96**

For illustrative purposes only.

To perform the assay the instructions for use provided with the kit have to be used.

Distributed by: **IBL International GmbH**
Flughafenstrasse 52a
22335 Hamburg, Germany

Phone: +49 (0)40-53 28 91-0 IBL@tecan.com
Fax: +49 (0)40-53 28 91-11 www.tecan.com/ibl

Always there for you



Code No. JP27776

Human APP β CTF Assay Kit - IBL**INTRODUCTION**

Alzheimer's disease (AD) was first reported by A. Alzheimer, a German neuropathologist in 1907 and is considered as a major factor of senile dementia. It is known that Amyloid β ($A\beta$; which is a major constituent of senile plaque) is cleaved from Amyloid Precursor Protein (APP; which exists in three main isoforms, APP695, APP751, and APP770) by β -secretase and subsequent γ -secretase (ref. 1). On the one hand, it has been also reported there is β CTF (β -secretase C-terminal fragment), fragments which are cleaved by β -secretase but not by γ -secretase.

It is thought that measuring β CTF, in addition to $A\beta$, soluble APP α and APP β , also has significance in research of the metabolic pathway of APP.

This ELISA can measure human APP β CTF in samples such as cultured cell lysate.

PRINCIPLE

This kit is a solid phase sandwich ELISA using 2 kinds of high specific antibodies. Tetra Methyl Benzidine (TMB) is used as a coloring agent (Chromogen). The strength of coloring is proportional to the quantities of Human APP β CTF.

MEASUREMENT RANGE

0.19 - 12 pmol/L

INTENDED USE

This IBL's assay kit is capable for the measurement of Human APP β CTF in cell lysate.

Preparation of the cultured cell lysate

1. Add IBLysis-I (IBL Lysate buffer #19022) to the cells pellet (2×10^7 indication).
2. Mix thoroughly with a mixer like Vortex.
3. Solubilize by rotation at 2-8°C for 30 minutes.
4. Centrifuge at 10,000 rpm for 10 minutes at 2-8°C.
5. Apply the supernatant diluted by "4, EIA buffer" as necessary.

KIT COMPONENT

1	Precoated plate : Anti-APP-C Rabbit IgG Affinity Purify	96Well x 1
2	Labeled antibody Conc.	
	: (30X) HRP conjugated Anti- Human $A\beta$ (N) (82E1) Mouse IgG MoAb Affinity Purify	0.4mL x 1
3	Standard : Synthetic peptide corresponding to Human APP β CTF	0.5mL x 2
4	EIA buffer*	30mL x 1
5	Solution for Labeled antibody*	12mL x 1
6	Chromogen : TMB solution	15mL x 1
7	Stop solution*	12mL x 1
8	Wash buffer Conc. *	50mL x 1

OPERATION MANUAL**1. Materials needed but not supplied**

- Plate reader (450nm)
- Graduated cylinder and beaker
- Refrigerator (as 4°C)
- Paper towel
- Washing bottle for precoated plate
- Disposable test tube for "2, Labeled antibody Conc." and "6, Chromogen"
- Micropipette and tip
- Deionized water
- Graph paper (log/log)
- Tube for dilution of Standard

2. Preparation

- 1) Preparation of wash buffer
"8, Wash buffer Conc." is a concentrated (40X) buffer. Adjust the temperature of "8, Washing buffer Conc." to room temperature and then, mix it gently and completely before use. Dilute 50 mL of "8, Wash buffer Conc." with 1,950 mL of deionized water and mix it. This is the wash buffer for use. This prepared wash buffer shall be stored in refrigerator and used within 2 weeks after dilution.

- 2) Preparation of Labeled antibody
"2, Labeled antibody Conc." is a concentrated (30X). Dilute "2, Labeled antibody Conc." with "5, Solution for Labeled antibody" in 30 times according to required quantity into a disposable test tube. Use this resulting solution as Labeled antibody.

Example)

In case you use one strip (8 well), the required quantity of Labeled antibody is 800 μ L. (Dilute 30 μ L of "2, Labeled antibody Conc." with 870 μ L of "5, Solution for Labeled antibody" and mix it. And use the resulting solution by 100 μ L in each well.)

This operation should be done just before the application of Labeled antibody.
The remaining "2, Labeled antibody Conc." should be stored at 4°C in firmly sealed vial.


- 3) Preparation of Standard
Put just 0.5 mL of deionized water into the vial of "3, Standard" and mix it gently and completely. This solution is 24 pmol/L human APP β CTF standard.
- 4) Dilution of Standard
Prepare 8 tubes for dilution of "3, Standard". Put 230 μ L each of "4, EIA buffer" into the tube.

Specify the following concentration of each tube."

Tube-1	12 pmol/L
Tube-2	6 pmol/L
Tube-3	3 pmol/L
Tube-4	1.5 pmol/L
Tube-5	0.75 pmol/L
Tube-6	0.38 pmol/L
Tube-7	0.19 pmol/L
Tube-8	0 pmol/L (Test Sample Blank)

Put 230 μ L of Standard solution into tube-1 and mix it gently. Then, put 230 μ L of tube-1 mixture into tube-2. Dilute two times standard solution in series to set up 7 points of diluted standard between 12 pmol/L and 0.19 pmol/L. Tube-8 is the test sample blank as 0 pmol/L.

See following picture.

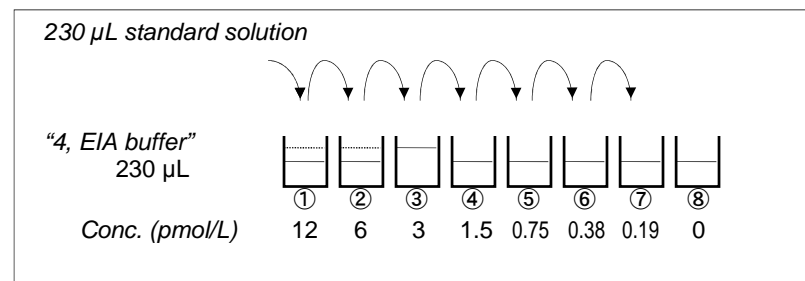

5) Dilution of test sample

Test sample should be diluted with "4, EIA buffer" as necessary.

If the quantity of human APP β CTF in samples may not be estimated in advance, the pre-assay with several different dilutions will be recommended to determine the proper dilution of samples.

3. Measurement procedure

All reagents shall be brought to room temperature approximately 30 minutes before use. Then mix it gently and completely before use. Make sure of no change in quality of the reagents. Standard curve shall be prepared simultaneously with the measurement of test samples.

Reagents	Test Sample	Standard	Test Sample Blank	Reagent Blank
	Test sample 100 μ L	Diluted standard (Tube 1-7) 100 μ L	EIA buffer (Tube-8) 100 μ L	EIA buffer 100 μ L
Incubation overnight at 4°C with plate lid				
4 times (wash buffer more than 350 μ L) *				
Labeled Antibody	100 μ L	100 μ L	100 μ L	-
Incubation for 60 minutes at 4°C with plate lid				
5 times (wash buffer more than 350 μ L) *				
Chromogen	100 μ L	100 μ L	100 μ L	100 μ L
Incubation for 30 minutes at room temperature (shielded)				
Stop solution	100 μ L	100 μ L	100 μ L	100 μ L
Read the plate at 450nm against a Reagent Blank within 30 minutes after addition of Stop solution.				

- 1) Determine wells for reagent blank. Put 100 μ L each of "4, EIA buffer" into the wells.
- 2) Determine wells for test sample blank, test sample and diluted standard. Then, put 100 μ L each of test sample blank (tube-8), test sample and dilutions of standard (tube-1-7) into the appropriate wells.
- 3) Incubate the precoated plate overnight at 4°C after covering it with plate lid.
- 4) Wash the plate with the prepared wash buffer and remove all liquid. *
- 5) Pipette 100 μ L of labeled antibody solution into the wells of test samples, diluted standard and test sample blank.
- 6) Incubate the precoated plate for 60 minutes at 4°C after covering it with plate lid.
- 7) Wash the plate with the prepared wash buffer and remove all liquid. *
- 8) Take the required quantity of "6, Chromogen" into a disposable test tube. Then, pipette 100 μ L from the test tube into the wells. Please do not return the rest of the test tube to "6, Chromogen" bottle to avoid contamination.
- 9) Incubate the precoated plate for 30 minutes at room temperature in the dark. The liquid will turn blue by addition of "6, Chromogen".
- 10) Pipette 100 μ L of "7, Stop solution" into the wells. Mix the liquid by tapping the side of precoated plate. The liquid will turn yellow by addition of "7, Stop solution".
- 11) Remove any dirt or drop of water on the bottom of the precoated plate and confirm there is no bubble on the surface of the liquid. Then, run the plate reader and conduct measurement at 450 nm against a reagent blank. The measurement shall be done within 30 minutes after addition of "7, Stop solution".

SPECIAL ATTENTION


- 1) Test samples should be measured soon after collection. For the storage of test samples, store them frozen and do not repeat freeze/thaw cycles. Thaw the test samples at a low temperature and mix them completely before measurement.
- 2) Test samples should be diluted with "4, EIA buffer", if the need arises.
- 3) Duplicate measurement of test samples and standard is recommended.
- 4) Use test samples in neutral pH range. The contaminations of organic solvent may affect the measurement.
- 5) Use only wash buffer contained in this kit for washing the precoated plate. Insufficient washing may lead to the failure in measurement.
- 6) Remove the wash buffer completely by tapping the precoated plate on paper towel. Do not wipe wells with paper towel.
- 7) "6, Chromogen" should be stored in the dark due to its sensitivity against light. "6, Chromogen" should be avoided contact with metals.
- 8) Measurement should be done within 30 minutes after addition of "7, Stop solution".

CALCULATION OF TEST RESULT

Subtract the absorbance of test sample blank from all data, including standards and unknown samples before plotting. Plot the subtracted absorbance of the standards against the standard concentration on log-log graph paper. Draw the best smooth curve through these points to construct the standard curve. Read the concentration for unknown samples from the standard curve.

Example of standard curve

Conc. (pmol/L)	Absorbance (450nm)
12	2.379
6	1.366
3	0.763
1.5	0.427
0.75	0.246
0.38	0.149
0.19	0.100
0 (Test Sample Blank)	0.047

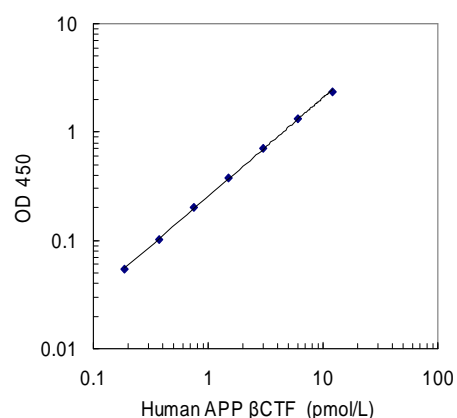

* The typical standard curve is shown above. This curve can not be used to derive test results. Please run a standard curve for each assay.

PERFORMANCE CHARACTERISTICS

1. Titer Assay (Samples with standard added are used.)

Specimen	Titer (X)	Measurement Value (pmol/L)	Theoretical Value (pmol/L)	%
NB-1 Cell lysate (IBLysis-I)	2	3.25	3.77	86.2
	4	1.63	1.94	84.0
	8	0.89	0.99	89.9
	16	0.51	0.47	108.5
Neuro2a Cell lysate (IBLysis-I)	4	1.23	1.58	77.8
	8	0.70	0.80	87.5
	16	0.36	0.41	87.8

IBLysis-I: Lysate buffer by IBL (# 19022)

2. Added Recovery Assay

Specimen	Theoretical Value (pmol/L)	Measurement Value (pmol/L)	%
IBLysis-I (x2)	3.00	2.92	97.3
	1.50	1.41	94.0
	0.75	0.71	94.7

3. Intra - Assay

Measurement Value (pmol/L)	SD value	CV value (%)	n
5.68	0.11	1.9	24
1.38	0.06	4.3	24
0.42	0.02	4.8	24

4. Inter - Assay

Measurement Value (pmol/L)	SD value	CV value (%)	n
5.39	0.23	4.3	5
1.32	0.05	3.8	5
0.44	0.01	2.3	5

5. Specificity


Compound	Cross Reactivity
Human APP βCTF	100%
Human sAPPβ-Wild Type	≤0.1%
Human sAPPβ-Swedish Type	≤0.1%
Human sAPPα	≤0.1%

6. Sensitivity

0.02 pmol/L

The sensitivity for this kit was determined using the guidelines under the National Committee for Clinical Laboratory Standards (NCCLS) Evaluation Protocols. (National Committee for Clinical Laboratory Standards Evaluation Protocols, SC1, (1989) Villanova, PA: NCCLS.)

7. Example of an experiment

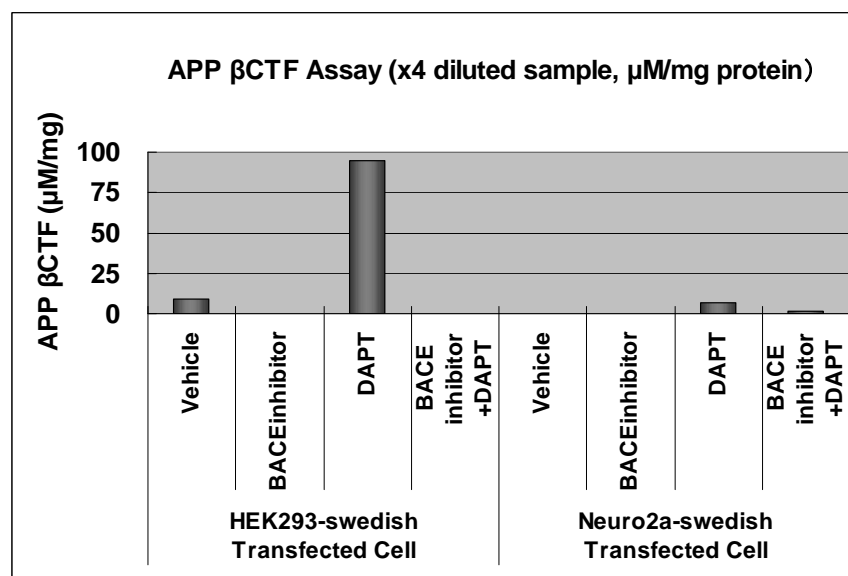

Data provided by Dr. Nobumasa Takasugi, Department of Neuropathology and Neuroscience, Graduate School of Pharmaceutical Sciences, University of Tokyo, Tokyo, Japan

PRECAUTION FOR INTENDED USE AND/OR HANDLING

- All reagents should be stored at 2 - 8°C. All reagents shall be brought to room temperature approximately 30 minutes before use.
- "3, Standard" is lyophilized products. Be careful to open this vial.
- "7, Stop solution" is a strong acid substance. Therefore, be careful not to have your skin and clothes contact "7, Stop solution" and pay attention to the disposal of "7, Stop solution".
- Dispose used materials after rinsing them with large quantity of water.
- Precipitation may occur in "2, Labeled antibody Conc.", however, there is no problem in the performance.
- Wash hands after handling reagents.
- Do not mix the reagents with the reagents from a different lot or kit.
- Do not use expired reagents.
- This kit is for research purpose only. Do not use for clinical diagnosis.

STORAGE AND THE TERM OF VALIDITY

Storage Condition: 2 - 8°C
The expiry date is specified on outer box.

REFERENCE

- Selkoe DJ. Normal and abnormal biology of the beta-amyloid precursor protein. Annu Rev Neurosci. 1994;17:489-517.

Version 2.












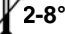



October 2016*

Amyloid β and APP related products:

Code No.	Product Name	Volume
27711	Human Amyloidβ (1-42) Assay Kit - IBL	96 Well
27712	Human Amyloidβ (1-42) (N) Assay Kit - IBL	96 Well
27713	Human Amyloidβ (1-40) Assay Kit - IBL	96 Well
27714	Human Amyloidβ (1-40) (N) Assay Kit - IBL	96 Well
27718	Human Amyloidβ (1-40) (FL) Assay Kit - IBL	96 Well
27720	Mouse/Rat Amyloidβ (1-40) High Specific Assay Kit - IBL	96 Well
27721*	Mouse/Rat Amyloidβ (1-42) Assay Kit - IBL*	96 Well
27716	Human Amyloidβ (N3pE-42) Assay Kit - IBL	96 Well
27418	Human Amyloidβ (N3pE-40) Assay Kit - IBL	96 Well
27729	Human Amyloidβ (1- x) Assay Kit - IBL	96 Well
27734*	Human sAPPα (highly sensitive) Assay Kit - IBL*	96 Well
27419*	Mouse/Rat sAPPα (highly sensitive) Assay Kit - IBL*	96 Well
27732*	Human sAPPβ-w (highly sensitive) Assay Kit - IBL*	96 Well
27733*	Human sAPPβ-sw (highly sensitive) Assay Kit - IBL*	96 Well
27776	Human APP βCTF Assay Kit - IBL	96 Well
19022	IBLysis-I (Lysate buffer)	50 mL 50x0.5mL

Made in Japan.

Symbols / Symbole / Symboles / Símbolos / Simboli / Símbolos / Σύμβολα

	Cat.-No.: / Kat.-Nr.: / No.- Cat.: / Cat.-No.: / N.-Cat.: / N.º Cat.: / Αριθμός-Κατ.:
	Lot-No.: / Chargen-Bez.: / No. Lot: / Lot-No.: / Lotto n.: / Lote N.º: / Αριθμός -Παραγωγή:
	Use by: / Verwendbar bis: / Utiliser à: / Usado por: / Da utilizzare entro:/ Usar até: / Χρησιμοποιείται από:
	No. of Tests: / Kitgröße: / Nb. de Tests: / No. de Determ.: / Quantità dei tests: / N.º de Testes: / Αριθμός εξετάσεων:
	Concentrate / Konzentrat / Concentré / Concentrar / Concentrato / Concentrado / Συμπύκνωμα
	Lyophilized / Lyophilisat / Lyophilisé / Liofilizado / Liofilizzato / Liofilizado / Λυοφιλοσμένο
	In Vitro Diagnostic Medical Device / In-vitro-Diagnostikum / Appareil Médical pour Diagnostics In Vitro / Dispositivo Médico para Diagnóstico In Vitro / Dispositivo Medico Diagnostico In vitro / Equipamento Médico de Diagnóstico In Vitro / Ιατρική συσκευή για In-Vitro Διάγνωση
	Contains biological material of human origin / Enthält biologisches Material menschlichen Ursprungs / Contient une substance biologique d'origine humaine / Contiene material biológico de origen humano / Contiene materiale biologico di origine umana / Contém material biológico de origem humana / Περιέχει βιολογικό υλικό ανθρώπινης προέλευσης
	Contains biological material of animal origin / Enthält biologisches Material tierischen Ursprungs / Contient une substance biologique d'origine animale / Contiene material biológico de origen animal / Contiene materiale biologico di origine animale / Contém material biológico de origem animal / Περιέχει βιολογικό υλικό ζωικής προέλευσης
	Unique Device Identification / Eindeutige Geräteerkennung / Identifiant de dispositif unique / Identificación única de producto / Identificatore univoco del dispositivo / Identificador de dispositivo único / Μοναδικός αναγνωριστικός κωδικός προϊόντος
	Read instructions before use / Arbeitsanleitung lesen / Lire la fiche technique avant emploi / Lea las instrucciones antes de usar / Leggere le istruzioni prima dell'uso / Ler as instruções antes de usar / Διαβάστε τις οδηγίες πριν την χρήση
	Keep away from heat or direct sun light / Vor Hitze und direkter Sonneneinstrahlung schützen / Garder à l'abri de la chaleur et de toute exposition lumineuse / Manténgase alejado del calor o la luz solar directa / Non esporre ai raggi solari / Manter longe do calor ou luz solar directa / Να φυλάσσεται μακριά από θερμότητα και άμεση επαφή με το φως του ηλίου
	Store at: / Lagern bei: / Stocker à: / Almacene a: / Armazena a: / Conservare a: / Armazena em: / Αποθήκευση στους:
	Store at: 2 - 8°C / Lagern bei: 2 - 8°C / Stocker à: 2 - 8°C / Almacene a: 2 - 8°C / Armazena a: 2 - 8°C / Conservare a: 2-8°C / Armazena em: 2-8°C / Αποθήκευση στους: 2-8°C
	Manufacturer: / Hersteller: / Fabricant: / Productor: / Fabricante: / Fabbricante: / Παραγωγός:
	Distributor: / Distributor: / Distributeur: / Distributor: / Distributore: / Distribuidor: / Διανομέας:
	Caution! / Vorsicht! / Attention! / ¡Precaución! / Attenzione! / Cuidado! / Προσοχή!
	Symbols of the kit components see MATERIALS SUPPLIED. Die Symbole der Komponenten sind im Kapitel KOMPONENTEN DES KITS beschrieben. Voir MATERIEL FOURNI pour les symboles des composants du kit. Símbolos de los componentes del juego de reactivos, vea MATERIALES SUMINISTRADOS. Per i simboli dei componenti del kit si veda COMPONENTI DEL KIT. Para símbolos dos componentes do kit ver MATERIAIS FORNECIDOS. Για τα σύμβολα των συστατικών του κιτ συμβουλευτείτε το ΠΑΡΕΧΟΜΕΝΑ ΥΛΙΚΑ.

Generic table, not all symbols are present in the product

COMPLAINTS: Complaints may be submitted initially written or vocal. Subsequently they need to be filed including the test performance and results in writing in case of analytical reasons.

WARRANTY: The product is warranted to be free from material defects within the specific shelf life and to comply with product specifications delivered with the product. The product must be used according to the Intended use, all instructions given in the instructions for use and within the product specific shelf life. Any modification of the test procedure or exchange or mixing of components of different lots could negatively affect the results. These cases invalidate any claim for replacement.

LIMITATION OF LIABILITY: IN ALL CIRCUMSTANCES THE EXTENT OF MANUFACTURER'S LIABILITY IS LIMITED TO THE PURCHASE PRICE OF THE KIT(S) IN QUESTION. IN NO EVENT SHALL MANUFACTURER BE LIABLE FOR ANY INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING DAMAGES FOR LOST PROFITS, LOST SALES, INJURY TO PERSON OR PROPERTY OR ANY OTHER INCIDENTAL OR CONSEQUENTIAL LOSS.

The labelling of hazardous substances is according to European directive.

For further country-specific classifications, please refer to the corresponding safety data sheet.



IBL International GmbH

Flughafenstrasse 52a
22335 Hamburg, Germany

Phone: +49 (0)40-53 28 91-0
Fax: +49 (0)40-53 28 91-11

IBL@tecan.com
www.tecan.com/ibl

Always there for you